

Dispenser 6 Magazine

Cat. 7158

Antibiotic discs dispenser with 6 cartridges

Practical information

Industry: Antimicrobial susceptibility testing

Principles and uses

The dispenser is a device for the manual application of antibiotic discs on the agar surface in Petri dish. It can only be used with the cartridges containing Laboratorios Conda's antibiotic discs.

Instructions for use

- Adjust the circular base of the dispenser according to the thickness of the agar.
- Insert the cartridge with antibiotic discs into the dispenser and make sure they are positioned correctly.
- Place the dispenser on a 90 mm Petri dish.
- Press the lever to the end and release. At this point, discs are placed on the agar medium.
- Dispenser must always perform a full cycle before it can work again.
- Note: If resistance is noticed when pressing the lever, stop immediately and make sure that there is no empty cartridge or that discs are not blocked.
- The dispenser locks automatically after running out of the discs in a cartridge. You must remove the empty cartridge to continue.

CLEANING

The dispenser should be cleaned with warm distilled water (you can add a few drops of liquid soap). If some parts of the dispenser are contaminated, they must be disinfected by immersing the dispenser in ethanol diluted in water (70% v / v) for 5 minutes, then washed with clean distilled water. Dry completely before using the dispenser again.

Do not use solvents such as acetone, chloroform, xylene and phenolic compounds or solutions of detergents.